

FD Rapid GolgiStain™ Kit

A complete Golgi-Cox staining system for the study of the morphology of neurons and glia

User Manual

PK 401/401A, Version 2019-01

FOR *IN VITRO* RESEARCH USE ONLY
not for diagnostic or other uses



FD NeuroTechnologies
Consulting & Services, Inc.

© 2002-2019 All Rights Reserved

Contents

I.	Introduction	4
II.	Kit Contents	5
III.	Materials Required but Not Included	5
IV.	Safety and Handling Precautions	6
V.	Tissue Preparation	6
VI.	Staining Procedure	13
VII.	Literature Cited	16
VIII.	Recent References Using This Kit	17

I. Introduction

The Golgi-Cox method (1987) has been one of the most effective techniques for studying the morphology of neuronal dendrites and dendritic spines. However, the reliability and time-consuming process of Golgi-Cox staining have been major obstacles to the widespread application of this technique. To overcome these shortcomings and to promote this invaluable technique, we have developed the FD Rapid GolgiStain™ Kit based on the principle of the methods described by Ramón-Moliner (1970) and Glaser and Van der Loos (1981).

This kit is reliable for visualizing morphological details of neurons, especially dendrites and dendritic spines. The kit is suitable for the analysis of various parameters of dendritic morphology—such as dendritic length and branching pattern and dendritic spine number, shape, and size—in different areas of both animal and postmortem human brains (Beggs, et al., 2003; An et. al., 2008; Milatovic et al., 2010; Koyama, 2013; Koyama et. al., 2015; Bicanic et al., 2017; Gstrein et al., 2018; for more references using this kit, visit www.fdneurotech.com). The kit can be used for processing large or small quantities of brain samples or sections all at once or over a period of time. A 40-min instructional video for tissue freezing, cryosectioning, and staining is provided (see www.fdneurotech.com).

The FD Rapid GolgiStain™ Kit should be used with fresh or shortly fixed brain tissue. Using with previously formalin-fixed or fresh-frozen brains is not recommended since satisfactory results with these types of samples are not typically obtained. Additionally, the kit does not work on any tissue that has already been cut, either free-floating sections or sections mounted on microscope slides (for more information, see Du, 2019, *Current Protocols in Neuroscience*, 88, e69. doi: 10.1002/cpns.69).

II. Kit Contents

Store at room temperature

	<u>PK401A</u>	<u>PK401</u>
Solution A	125 ml	250 ml
Solution B	125 ml	250 ml
Solution C	125 ml x 2	250 ml x 2
Solution D	125 ml	250 ml
Solution E	125 ml	250 ml
Glass specimen retriever	2	2
Natural hair paintbrush	2	2
Dropping bottle	1	1
Plastic forceps	1	1
User manual	1	1

III. Materials Required but Not Included

1. Double distilled or Milli-Q water
2. Plastic/glass tubes or vials
3. Histological supplies and equipment:
 - Gelatin-coated microscope slides (Cat. #PO101)
 - Coverslips
 - Staining jars
 - Ethanol
 - Xylene
 - Eukitt[®] Quick-Hardening Mounting medium
 - Cryostat or vibratome
 - Upright bright-field microscope

Eukitt is a registered trademark of O. Kindler GmbH & Co.

IV. Safety and Handling Precautions*

1. FD Rapid GolgiStain™ Kit is made for *in vitro* research use only and not for drug, diagnostic or other uses.
2. The kit contains reagents that are toxic and harmful in contact with skin or by inhalation and may be fatal if ingested. Do not pipette by mouth. Avoid inhalation and contact with skin and eyes. In case of contact, wash immediately with generous amounts of water and seek medical advice. If swallowed, wash out mouth with water and immediately call a physician.
3. Perform experiment under a chemical hood. **Wear suitable protective clothing, gloves and eye/face protection while handling kit reagents.** Wash hands thoroughly after performing the experiment.

*Material safety data sheet (MSDS) is available at www.fdtypeurotech.com.

V. Tissue Preparation

Please read the following instructions before using this kit.

*See also an instructional video for tissue freezing and cryosectioning at www.fdtypeurotech.com or *Current Protocols in Neuroscience*, 88, e69. doi: 10.1002/cpns.69.*

- All containers (plastic preferred) to be used should be cleansed and rinsed with distilled water.
- Do not use metal implements whenever Solutions A and B are present.
- Keep containers tightly closed at all times.

- Tissues, including sections treated with Solutions A and B, should be protected from light whenever possible.
 - The following procedure should be performed at room temperature unless specifically indicated.
1. Prepare impregnation solution (Solution A/B) by mixing equal volumes of Solutions A and B at least 24 hours prior to use and leave unstirred.

The impregnation solution should be stored at room temperature in the dark before use and used within 4 weeks.

CAUTION: Solutions A and B (which contain mercuric chloride, potassium dichromate, and potassium chromate) are toxic if they come into contact with skin and may be fatal if swallowed. Do not breathe vapor or fumes from these solutions.

CAUTION: Do not pour the waste of Solutions A and B into the sink. Collect waste in a waste container, and contact the local safety office or a licensed professional waste disposal service to dispose of this material.

2. Deeply anesthetize experimental animal with IACUC-approved methods before sacrificing.

Do not perfuse animals unless absolutely necessary. If perfusion is necessary, tissue must NOT be postfixed (see step 3 for more details).

3. *Optional:* Perfuse animal with 4% paraformaldehyde in 0.1 M phosphate buffer (pH 7.4) according to the following suggested protocol (e.g., for both Golgi-Cox impregnation and immunostaining).

- a. Perfuse animal with 4% paraformaldehyde in 0.1 M phosphate buffer (pH 7.4) for 4 min.
 - b. Remove brain from skull immediately after perfusion.
 - c. Divide brain (e.g., mouse or rat brains) into 2 hemispheres or 5-mm thick blocks.
 - d. For Golgi-Cox staining, continue to step 5.
DO NOT POSTFIX TISSUE!
 - e. For immunostaining, continue to process tissue according to desired protocol.
4. Remove animal brain (or postmortem human brain tissue) from the skull as quickly as possible. Be sure to handle carefully to avoid damaging or pressing the tissue.

Large brain specimens, including rat brains, should be divided with a sharp blade into blocks of approximately 5 to 10 mm thickness. For example, a rat brain may be divided either sagittally into 2 hemispheres or coronally into 2 to 3 blocks depending on the regions of interest.

5. Rinse tissue quickly in double distilled or Milli-Q water to remove blood from the surface.
6. Immerse tissue in impregnation solution (Solution A/B; from step 1), and store at room temperature for 2 weeks in the dark. Replace impregnation solution after the first 6 hr of immersion or on the next day. Gently swirl (do not shake!) the tissue container side to side for a few seconds at least twice a week during the impregnation period.

Use at least 5 ml impregnation solution for each cm³ of tissue. However, it is strongly recommended to use 1.5

times the minimum amount of impregnation solution (e.g., for a mouse brain, instead of using 5 ml, use 7.5 ml). Note that using a lower volume of impregnation solution may decrease the intensity and reliability of staining. The impregnation solution should be prepared at least 24 hours prior to use and left unstirred. It is important to use the top part of the solution that is free of precipitate.

In most cases, 2 weeks of impregnation is satisfactory. However, variations in type and size of tissue may require shorter or longer impregnation to obtain the best results. The optimal time should be obtained by trial for each type of tissue, but 3 weeks should be sufficient for most tissues. Note that prolonging the impregnation may increase background staining.

7. Transfer tissue into Solution C, and store at room temperature in the dark for at least 72 hr (up to 1 week). Replace Solution C at least once after the first 24 hours of immersion or on the next day.

The amount of Solution C should be at least 7 times the tissue volume. Using a lower volume of Solution C may result in higher background staining and nonspecific precipitate.

8. Freeze tissue for sectioning with a cryostat or sliding microtome. For example, to rapidly freeze tissue:
 - a. Place tissue in a plastic spoon, and slowly dip into isopentane that has been precooled with dry ice.

The temperature should be kept below -70°C , and the dipping should take about 1 min for a mouse or rat brain (the slower the better).

- b. After the tissue is completely immersed in isopentane, place on dry ice for another minute to ensure that the tissue is well frozen.

Do not let the tissue thaw before the sections are cut.

To prevent damage to tissue caused by ice crystal formation and to preserve the best possible cell morphology, the tissue must be properly frozen before sectioning.

9. Cut 100- to 200- μm sections using a cryostat (recommended) at -20°C to -23°C .

If the cryostat has only one temperature control, set the cryostat temperature to -23°C at least 4 hrs before cutting. If the cryostat has 2 temperature settings, set the chamber temperature 1°C colder than that of the specimen head. In most cases, -23°C is satisfactory; however, variations in type of cryostat and tissue may require a higher or lower chamber temperature in order to cut sections smoothly and without shattering.

- a. Mount tissue on the specimen disc with distilled water, but make sure that the tissue does not thaw (may be done on dry ice).

Tissue may also be mounted with any type of tissue freezing medium, including optimal cutting temperature (OCT) compound, but avoid cutting through the freezing medium. Do not embed the tissue in OCT. If the tissue must be embedded for cutting, use TFM™ tissue freezing medium.

To help prevent sections from sticking to the blade, the brain surface may be coated with a thin layer of distilled water (ice) using a paintbrush. Make sure that the brain

does not thaw at any time. The brain can be coated either before or after mounting on the specimen disc. When coating with water, the brain should be very cold so that the water freezes immediately. To coat the brain before mounting on specimen disc, first place the brain directly on dry ice. Then, dip a paintbrush in cold, deionized water, and brush the surface of the brain to apply a thin layer of water using brush strokes that are fast enough to avoid paint brush hairs sticking to the brain.

- b. After tissue is mounted on the specimen disc, place tissue on dry ice for 10 min so that both the tissue and specimen disc touch the dry ice.
- c. Immediately install specimen disc with tissue on the cryostat, and wait 3 to 5 min before cutting 1 or 2 sections.

Do not use the anti-roll plate.

- d. Set cryostat to the thickness of sections to be cut, and cut sections.

Each type of cryostat may vary, but all types should be able to cut thick sections. If the cryostat does not have the option to set the section thickness to $>60\ \mu\text{m}$, such as for 100- μm sections (or thicker), cut sections as follows: (1) Set cryostat to 50- μm section thickness. (2) Roll cryostat counterclockwise slowly, and stop when the tissue almost touches the blade (or knife). (3) Roll cryostat clockwise to cut the tissue. If the tissue is too cold (e.g., sections show cracks that are parallel to the blade), wait 1 to 2 min before the next trial. Repeat this step until high-quality sections are achieved, and then continue cutting. (4) Repeat to cut each section.

Optional: Cut 100- to 200- μm sections with other types of microtomes, such as vibratome, sliding microtome, or rotary microtome.

If using a vibratome, the impregnated brain should be embedded in agarose or gelatin. However, for collecting sections, the cutting chamber must be filled with Solution C, otherwise sections may crack upon drying.

If using a sliding microtome, both the stage and blade need to be maintained at low temperature ($<0^{\circ}\text{C}$).

If using a rotary microtome, tissue must be embedded in paraffin after immersion in Solution C (step 7).

10. Mount sections on gelatin-coated microscope slides using Solution C.

- a. Apply a few drops of Solution C (dropper bottle provided) onto a gelatin-coated microscope slide.
- b. Transfer sections with a glass specimen retriever (provided) onto the slide.
- c. After moving each section to an ideal location with a paintbrush, remove excess Solution C from the slide using a Pasteur pipette and a strip of filter paper if necessary.

Any solution on the slide must be removed as much as possible; otherwise, sections may fall off the slides during staining.

- d. Allow sections to dry naturally overnight at room temperature. Do not use a fan or hot plate.

Sections should be stained as soon as possible after drying

overnight, but unstained sections may be stored in a slide box at room temperature in the dark for up to 3 days before staining.

To prevent sections from falling off slides, properly prepared gelatin-coated microscope slides must be used (e.g., FD NeuroTechnologies, Cat. No. PO101).

VI. Staining Procedure

Do not let sections dry out between any steps during the staining or while coverslipping!

*See also an instructional video for staining at www.fdnurotech.com or *Current Protocols in Neuroscience*, 88, e69. doi: 10.1002/cpns.69.*

1. Prepare staining solution (Solution D/E), which consists of 1 part Solution D, 1 part Solution E, and 2 parts double distilled or Milli-Q water. For example, mix the following solutions in the order listed: 10 ml Solution D, 10 ml Solution E, and 20 ml double distilled water.
 - *The staining solution should be prepared immediately before use and may be used for up to 100 sections (e.g., mouse brain) per 100 ml, depending on the size of sections.*
 - *Keep the bottle and staining jar containing the staining solution covered to prevent evaporation.*
2. Rinse sections in double distilled or Milli-Q water 2 times for 4 min each rinse. Replace distilled water after each use.

To help prevent sections from falling off slides, cold double distilled or Milli-Q water that has been stored at 4°C may be used.

3. Place sections in the staining solution (Solution D/E) for 10 min.

Sections should be stained in a glass or plastic staining jar with sufficient solution. Do not stain sections on microscope slides.

The staining solution should be stirred frequently during staining.

4. Rinse sections in double distilled or Milli-Q water 2 times for 4 min each rinse. Replace distilled water after each use.

To help prevent sections from falling off slides, cold double distilled or Milli-Q water that has been stored at 4°C may be used.

5. Optional: Counterstain Golgi impregnated sections with cresyl violet.

***Note:** Counterstaining allows not only determination of the precise localization of impregnated neurons but also visualization of anatomic landmarks that would not be evident in a non-counterstained section. However, some fine dendritic branches and dendritic spines of impregnated neurons may be obscured by cresyl violet-stained cellular elements.*

Do not let sections dry out between any steps described below.

a. Continue to rinse sections in double distilled or Milli-Q water 2 times for 4 min each rinse. Replace distilled water after each use.

b. Place sections in 50% ethanol for 5 min.

c. Stain sections in cresyl violet solution for 20 to 30 min.

Staining time may be increased or decreased depending on the desired intensity and the concentration of cresyl violet solution used.

d. Dip sections in double distilled or Milli-Q water 3 times. Replace distilled water after each use.

e. Dehydrate sections in sequential rinses of 50%, 75%, and 95% ethanol, 30 to 50 sec each rinse, depending on the desired intensity. Do not skip any step.

The staining intensity of cellular elements and background decreases quickly in these solutions.

f. Continue to step 7.

6. Dehydrate sections in sequential rinses of 50%, 75%, and 95% ethanol, 4 min each rinse. Do not skip any step.

7. Dehydrate sections in 100% ethanol 4 times for 4 min each rinse.

Incubation may be prolonged if the section thickness is >150 μ m.

8. Clear sections in xylene 3 times for 4 min each rinse. *Sections may be temporarily stored in xylene for a few hours before overslipping.*

9. Coverslip sections with Eukitt[®] Quick-hardening mounting medium or Permout[®] mounting medium if the former is not available.

Our recent tests indicate that the Eukitt[®] Quick-hardening mounting medium minimizes the development of nonspecific spherical-shaped precipitates in the background after prolonged storage and is therefore recommended for use with the FD Rapid GolgiStain[™] Kit.

10. Golgi-stained sections should be stored at room temperature and protected from light.

Golgi-stained sections should be imaged or analyzed as soon as possible. Generally, sections can be stored at room temperature in the dark for 1 year without decrease in staining intensity.

VII. Literature Cited

- An, J. J., Gharami, K., Liao, G., Woo, N. H., Lau, A. G., Vanevski, F., Torre, E. R., Jones, K. R., Feng, Y., Lu, B., Xu, B. (2008). Distinct Role of Long 3' UTR BDNF mRNA in Spine Morphology and Synaptic Plasticity in Hippocampal Neurons. *Cell*, 134(1), 175-187. doi:10.1016/j.cell.2008.05.045
- Beggs, H. E., Schahin-Reed, D., Zang, K., Goebbels, S., Nave, K., Gorski, J., Jones, K. R., Sretavan, D., Reichardt, L. F. (2003). FAK Deficiency in Cells Contributing to the Basal Lamina Results in Cortical Abnormalities Resembling Congenital Muscular Dystrophies. *Neuron*, 40(3), 501-514. doi:10.1016/s0896-6273(03)00666-4
- Bicanic, I., Hladnik, A., & Petanjek, Z. (2017). A Quantitative Golgi Study of Dendritic Morphology in the Mice Striatum Medium Spiny Neurons. *Frontiers in Neuroanatomy*, 11. doi:10.3389/fnana.2017.00037
- Corsi, P. (1987). Camillo Golgi's morphological approach to neuroanatomy. In Masland R. L., Portera-Sanchez A. and Toffano G. (Ed.), *Neuroplasticity: A new therapeutic tool in the CNS pathology* (pp.1-7). Padova: Liviviana Press.
- Glaser, E. M., & Van der Loos, H. (1981). Analysis of thick brain sections by obverse—reverse computer microscopy: Application of a new, high clarity Golgi—Nissl stain. *Journal of Neuroscience Methods*, 4(2), 117-125. doi:10.1016/0165-0270(81)90045-5

- Gstrein, T., Edwards, A., Pristoupilová, A., Leca, I., Breuss, M., Pilat-Carotta, S., Hansen, A. H., Tripathy, R., Traunbauer, A. K., Hochstoeger, T., Rosoklija, G., Repic, M., Landler, L., Stránecký, V., Dürmberger, G., Keane, T. M., Zuber, J., Adams, D. J., Flint, J., Honzik, T., Gut, M., Beltran, S., Mechtler, K., Sherr, E., Kmoch, S., Gut, I., Keays, D. A. (2018). Mutations in Vps15 perturb neuronal migration in mice and are associated with neurodevelopmental disease in humans. *Nature Neuroscience*, 21(2), 207-217. doi:10.1038/s41593-017-0053-5
- Koyama, Y. (2013). The unending fascination with the Golgi method. *OA Anatomy*, 1(3). doi:10.13172/2052-7829-1-3-848
- Koyama, Y., Hattori, T., Nishida, T., Hori, O., & Tohyama, M. (2015). Alterations in dendrite and spine morphology of cortical pyramidal neurons in DISC1-binding zinc finger protein (DBZ) knockout mice. *Frontiers in Neuroanatomy*, 9. doi:10.3389/fnana.2015.00052
- Milatovic, D., Montine, T. J., Zaja-Milatovic, S., Madison, J. L., Bowman, A. B., & Aschner, M. (2010). Morphometric Analysis in Neurodegenerative Disorders. *Current Protocols in Toxicology*. doi:10.1002/0471140856.tx1216s46
- Ramón-Moliner E. (1970). The Golgi-Cox technique. In Nauta W.J.H and Ebbsen S.O.E. (Ed.), *Contemporary research methods in neuroanatomy: With 190 figures* (pp.32-55). Berlin: Springer.

VIII. Recent References Using This Kit

- Ano, Y., Ohya, R., Kita, M., Taniguchi, Y. and Kondo, K., (2019). Theaflavins Improve Memory Impairment and Depression-Like Behavior by Regulating Microglial Activation. *Molecules*, 24(3), p.467.
- Apaijai, N., Moisesu, D.M., Palee, S., McSweeney, C.M., Saiyasit, N., Maneechote, C., Boonnag, C., Chattipakorn, N. and Chattipakorn, S.C., (2019). Pretreatment With PCSK9 Inhibitor Protects the Brain Against Cardiac Ischemia/Reperfusion Injury Through a Reduction of Neuronal Inflammation and Amyloid Beta Aggregation. *Journal of the American Heart Association*, 8(2), p.e010838.
- Beckelman, B.C., Yang, W., Kasica, N.P., Zimmermann, H.R., Zhou, X., Keene, C.D., Ryazanov, A.G. and Ma, T., (2019). Genetic reduction of eEF2 kinase alleviates pathophysiology in Alzheimer's disease model mice. *The Journal of clinical investigation*, 129(2), 820-833.
- Bello-Medina, P.C., Prado-Alcalá, R.A. and Rivas-Arancibia, S., (2019). Effect of ozone exposure on dendritic spines of CA1 pyramidal neurons of the dorsal hippocampus and on object-place recognition memory in rats. *Neuroscience*. PMID:30685541
- Boros, B.D., Greathouse, K.M., Gearing, M. and Herskowitz, J.H., (2019). Dendritic spine remodeling accompanies Alzheimer's disease pathology and genetic susceptibility in cognitively normal aging. *Neurobiology of aging*, 73, pp.92-103.
- Bowman, R.E., Hagedorn, J., Madden, E. and Frankfurt, M., (2019). Effects of adolescent Bisphenol-A exposure on memory and spine density in ovariectomized female rats: Adolescence vs adulthood. *Hormones and behavior*, 107, pp.26-34.
- Chen, Y.F., Chen, Z.X., Wang, R.H., Shi, Y.W., Xue, L., Wang, X.G. and Zhao, H., (2019). Knockdown of CLC-3 in the hippocampal CA1 impairs contextual fear memory. *Progress in Neuro-Psychopharmacology and Biological Psychiatry*, 89, pp.132-145.
- Czechowska, N., van Rienen, A., Lang, F., Eiberger, B. and Baader, S.L., (2019). An update on the Golgi staining technique improving cerebellar cell type specificity. *Histochemistry and cell biology*, PMID:30607497
- Daemi, F., Aumann, T., Biemond, M., Wood, R.J., Hoffmann, E., Pasquet, J., Vessey, K.A., Nicholson, M., Fletcher, J.L., Gonsalvez, D.G. and Gunnarsen, J.M., (2019). Neuronal TrkB Drives Oligodendrocyte Production and Central Myelination. *Available at SSRN 3318928*.
- Dong, Q., Teng, S.W., Wang, Y., Qin, F., Li, Y., Ai, L.L. and Yu, H., (2019). Sitagliptin protects the cognition function of the Alzheimer's disease mice through activating glucagon-like peptide-1 and BDNF-TrkB signalings. *Neuroscience letters*, 696, pp.184-190.
- Hong, Y., Jin, Y., Park, K., Choi, J., Kang, H., Lee, S.R. and Hong, Y., (2019). Elevated Serum Melatonin under Constant Darkness Enhances Neural Repair in Spinal Cord Injury through Regulation of Circadian Clock Proteins Expression. *Journal of clinical medicine*, 8(2), p.135.

12. Hu, Y., Chen, W., Wu, L., Jiang, L., Liang, N., Tan, L., Liang, M. and Tang, N., (2019). TGF- β 1 Restores Hippocampal Synaptic Plasticity and Memory in Alzheimer Model via the PI3K/Akt/Wnt/ β -Catenin Signaling Pathway. *Journal of Molecular Neuroscience*, 67(1):142-149.
13. Kawase, H., Ago, Y., Naito, M., Higuchi, M., Hara, Y., Hasebe, S., Tsukada, S., Kasai, A., Nakazawa, T., Mishina, T. and Kouji, H., (2019). mS-11, a mimetic of the mSin3-binding helix in NR5F, ameliorates social interaction deficits in a prenatal valproic acid-induced autism mouse model. *Pharmacology Biochemistry and Behavior*, 176, pp.1-5.
14. Liu, B., Ou, G., Chen, Y. and Zhang, J., (2019). Inhibition of protein tyrosine phosphatase 1B protects against sevoflurane-induced neurotoxicity mediated by ER stress in developing brain. *Brain research bulletin*, 146, pp.28-39.
15. Noh, K., Lee, H., Choi, T.Y., Joo, Y., Kim, S.J., Kim, H., Kim, J.Y., Jahng, J.W., Lee, S., Choi, S.Y. and Lee, S.J., (2019). Negr1 controls adult hippocampal neurogenesis and affective behaviors. *Molecular Psychiatry*, PMID:30651602
16. Parato, J., Shen, H. and Smith, S.S., (2019). α 4 β 8 GABAA Receptors Trigger Synaptic Pruning and Reduce Dendritic Length of Female Mouse CA3 Hippocampal Pyramidal Cells at Puberty. *Neuroscience*, 398, pp.23-36.
17. Parra, A., Baker, C.A. and Bolton, M.M., (2019). Regional Specialization of Pyramidal Neuron Morphology and Physiology in the Tree Shrew Neocortex. *Cerebral Cortex*. PMID:30715235
18. Peng, J., Pang, J., Huang, L., Enkhjargal, B., Zhang, T., Mo, J., Wu, P., Xu, W., Zuo, Y., Peng, J. and Zuo, G., (2019). LRP1 activation attenuates white matter injury by modulating microglial polarization through Shc1/PI3K/Akt pathway after subarachnoid hemorrhage in rats. *Redox Biology*, p.101121.
19. Schaefer, M.L., Wang, M., Perez, P.J., Peralta, W.C., Xu, J. and Johns, R.A., (2019). Nitric Oxide Donor Prevents Neonatal Isoflurane-induced Impairments in Synaptic Plasticity and Memory. *Anesthesiology: The Journal of the American Society of Anesthesiologists*, 130(2), pp.247-262.
20. Sripetchwandee, J., Pintana, H., Sa-nguanmoo, P., Boonnag, C., Pratchayasakul, W., Chattipakorn, N. and Chattipakorn, S.C., (2019). Comparative effects of sex hormone deprivation on the brain of insulin-resistant rats. *Journal of Endocrinology*, PMID: 30689543.
21. Thelen, C., Flaherty, E., Saurine, J., Sens, J., Mohamed, S. and Pitychoutis, P.M., (2019). Sex Differences in the Temporal Neuromolecular and Synaptogenic Effects of the Rapid-acting Antidepressant Drug Ketamine in the Mouse Brain. *Neuroscience*, 398, pp.182-192.
22. Trzesniewski, J., Altmann, S., Jäger, L. and Kapfhammer, J.P., (2019). Reduced Purkinje cell size is compatible with near normal morphology and function of the cerebellar cortex in a mouse model of spinocerebellar ataxia. *Experimental neurology*, 311, pp.205-212.
23. Wang, L., Pang, K., Han, K., Adamski, C.J., Wang, W., He, L., Lai, J.K., Bondar, V.V., Duman, J.G., Richman, R. and Toliaas, K.F., (2019). An autism-linked missense mutation in SHANK3 reveals the modularity of Shank3 function. *Molecular psychiatry*, PMID:30610205
24. Wang, X.M., Pan, W., Xu, N., Zhou, Z.Q., Zhang, G.F. and Shen, J.C., (2019). Environmental enrichment improves long-term memory impairment and aberrant synaptic plasticity by BDNF/TrkB signaling in nerve-injured mice. *Neuroscience letters*, 694, pp.93-98.
25. Xie, Q., Cheng, J., Pan, G., Wu, S., Hu, Q., Jiang, H., Wang, Y., Xiong, J., Pang, Q. and Chen, X., (2019). Treadmill exercise ameliorates focal cerebral ischemia/reperfusion-induced neurological deficit by promoting dendritic modification and synaptic plasticity via upregulating caveolin-1/VEGF signaling pathways. *Experimental neurology*, 313, pp.60-78.
26. Yan, X.S., Yang, Z.J., Jia, J.X., Song, W., Fang, X., Cai, Z.P., Huo, D.S. and Wang, H., (2019). Protective mechanism of testosterone on cognitive impairment in a rat model of Alzheimer's disease. *Neural regeneration research*, 14(4), p.649-657.
27. Yu, M., Chen, X., Liu, J., Ma, Q., Zhuo, Z., Chen, H., Zhou, L., Yang, S., Zheng, L., Ning, C. and Xu, J., (2019). Gallic acid disruption of A β 1-42 aggregation rescues cognitive decline of APP/PS1 double transgenic mouse. *Neurobiology of disease*, 124, pp.67-80.
28. Yu, P., Venkat, P., Chopp, M., Zacharek, A., Shen, Y., Liang, L., Landschoot-Ward, J., Liu, Z., Jiang, R. and Chen, J., (2019). Deficiency of tPA Exacerbates White Matter Damage, Neuroinflammation, Glymphatic Dysfunction and Cognitive Dysfunction in Aging Mice. *Aging and disease*, 10.14336/AD.2018.0816.
29. Zhang, L., Dong, H., Si, Y., Wu, N., Cao, H., Mei, B. and Meng, B., (2019). miR-125b promotes tau phosphorylation by targeting the neural cell adhesion molecule in neuropathological progression. *Neurobiology of aging*, 73, pp.41-49.
30. Zhang, W.J., Cao, W.Y., Huang, Y.Q., Cui, Y.H., Tu, B.X., Wang, L.F., Zou, G.J., Liu, Y., Hu, Z.L., Hu, R. and Li, C.Q., (2019). The Role of miR-150 in Stress-Induced Anxiety-Like Behavior in Mice. *Neurotoxicity research*, 35(1), pp.160-172.
31. Zhu, R., Pan, Y.H., Sun, L., Zhang, T., Wang, C., Ye, S., Yang, N., Lu, T., Wisniewski, T., Tang, S. and Zhang, W., (2019). ADAMTS18 Deficiency Affects Neuronal Morphogenesis and Reduces the Levels of Depression-like Behaviors in Mice. *Neuroscience*, 399, pp.53-64.
32. Acharjee, S., Verbeek, M., Gomez, C.D., Bisht, K., Lee, B., Benoit, L., Sharkey, K.A., Benediktsson, A., Tremblay, M.E. and Pittman, Q.J., (2018). Reduced microglial activity and enhanced glutamate transmission in the basolateral amygdala in early CNS autoimmunity. *Journal of Neuroscience*, 38(42), pp.9019-9033.

33. Aguayo, F.I., Tejos-Bravo, M., Díaz-Véliz, G., Pacheco, A., García-Rojo, G., Corrales, W., Olave, F.A., Aliaga, E., Ulloa, J.L., Avalos, A.M. and Román-Albasini, L. (2018). Hippocampal Memory Recovery After Acute Stress: A Behavioral, Morphological and Molecular Study. *Frontiers in molecular neuroscience*, 11, 283.
34. Al-Absi, A.R., Christensen, H.S., Sanchez, C. and Nyengaard, J.R., (2018). Evaluation of semi-automatic 3D reconstruction for studying geometry of dendritic spines. *Journal of chemical neuroanatomy*, 94, pp.119-124.
35. Alvarez-Periel, E., Puigdemívol, M., Brito, V., Plattner, F., Bibb, J.A., Alberch, J. and Ginés, S. (2018). Cdk5 Contributes to Huntington's disease learning and memory deficits via modulation of brain region-specific substrates. *Molecular neurobiology*, 55(8):6250-6268.
36. Alves, M.J., Goksel, M., Kaya, B., Mostafa, H., Gygli, P., Stephens, J., Fair, S., Otero, J.J. and Czeisler, C.M. (2018). CCNA2 Ablation in Aged Mice Results in Abnormal rRNA Granule Accumulation in Hippocampus. *The American Journal of Pathology*, PubMed PMID: 30579783.
37. Arbogast, T., Razaz, P., Ellegood, J., McKinstry, S., Erdin, S., Currall, B., Anechych, T., Lerch, J.P., Qiu, L.R., Rodriguiz, R.M. and Mark Henkelman, R. (2018). Kctd13-deficient mice display short-term memory impairment and sex-dependent genetic interactions. *Human Molecular Genetics*, PubMed PMID: 30590535.
38. Ashokan, A., Lim, J.W.H., Hang, N. and Mitra, R. (2018). Complex housing causes a robust increase in dendritic complexity and spine density of medial prefrontal cortical neurons. *Scientific reports*, 8(1):7308.
39. Ayata, P., Badimon, A., Strasburger, H.J., Duff, M.K., Montgomery, S.E., Loh, Y.H.E., Ebert, A., Pimenova, A.A., Ramirez, B.R., Chan, A.T. and Sullivan, J.M., (2018). Epigenetic regulation of brain region-specific microglia clearance activity. *Nature neuroscience*, 21(8), p.1049.
40. Aytan, N., Choi, J.K., Carreras, I., Crabtree, L., Nguyen, B., Lehar, M., Blusztajn, J.K., Jenkins, B.G. and Dedeoglu, A. (2018). Protective effects of 7, 8-dihydroxyflavone on neuropathological and neurochemical changes in a mouse model of Alzheimer's disease. *European journal of pharmacology*, 828:9-17.
41. Banks, G., Lassi, G., Hoerder-Suabedissen, A., Tinarelli, F., Simon, M.M., Wilcox, A., Lau, P., Lawson, T.N., Johnson, S., Rutman, A. and Sweeting, M. (2018). A missense mutation in Katnall1 underlies behavioural, neurological and ciliary anomalies. *Molecular psychiatry*, 23(3):713.
42. Belanger, A.M., Przybylska, M., Geffras, E., Furgerson, M., Geller, S., Kloss, A., Cheng, S.H., Zhu, Y. and Yew, N.S. (2018). Inhibiting neutral amino acid transport for the treatment of phenylketonuria. *JCI insight*, 3 (14).
43. Blair, J.A. (2018). *Luteinizing hormone in the central nervous system: a direct role in learning and memory (Doctoral dissertation, Kent State University)*.
44. Blanco-Suarez, E., Liu, T.F., Kopelevich, A. and Allen, N.J. (2018). Astrocyte-Secreted Chordin-like 1 Drives Synapse Maturation and Limits Plasticity by Increasing Synaptic GluA2 AMPA Receptors. *Neuron*, 100 (5):1116-1132.
45. Bracke, A., 2018. Investigation of the effects of obesity on hippocampal morphometry and neuronal plasticity as well as behavior in a leptin-deficient mouse model (*Doctoral dissertation, Faculty of Mathematics and Natural Sciences of the Ernst Moritz Arndt University Greifswald, Germany*).
46. Bracke, A., Schäfer, S., und Halbach, V.V.B., Klempin, F., Bente, K., Bracke, K., Staar, D., van den Brandt, J., Harzsch, S., Bader, M. and Wenzel, U.O. (2018). ATP6AP2 over-expression causes morphological alterations in the hippocampus and in hippocampus-related behaviour. *Brain Structure and Function*, 223(5):2287-2302.
47. Carr, H., Alexander, T.C., Groves, T., Kiffer, F., Wang, J., Price, E., Boerma, M. and Allen, A.R. (2018). Early effects of 160 radiation on neuronal morphology and cognition in a murine model. *Life sciences in space research*, 17:63-73.
48. Chai, H., Liu, B., Zhan, H., Li, X., He, Z., Ye, J., Guo, Q., Chen, J., Zhang, J. and Li, S. (2018). Antidepressant Effects of Rhodomrytone in Mice with Chronic Unpredictable Mild Stress-Induced Depression. *International Journal of Neuropsychopharmacology*, PubMed PMID: 30407505.
49. Chatterjee, S., Cassel, R., Schneider-Anthony, A., Merienne, K., Cosquer, B., Tzeplaeff, L., Sinha, S.H., Kumar, M., Chaturbedy, P., Eswaremoorthy, M. and Le Gras, S. (2018). Reinstating plasticity and memory in a tauopathy mouse model with an acetyltransferase activator. *EMBO molecular medicine*, 10(11):e8587.
50. Chen, J., Niu, Q., Xia, T., Zhou, G., Li, P., Zhao, Q., Xu, C., Dong, L., Zhang, S. and Wang, A. (2018). ERK1/2-mediated disruption of BDNF-TrkB signaling causes synaptic impairment contributing to fluoride-induced developmental neurotoxicity. *Toxicology*, 410:222-230.
51. Chen, Y.W., Akad, A., Aderogba, R., Chowdhury, T.G. and Aoki, C. (2018). Dendrites of the dorsal and ventral hippocampal CA1 pyramidal neurons of singly housed female rats exhibit lamina-specific growths and retractions during adolescence that are responsive to pair housing. *Synapse*, 72(7):e22034.
52. Cheng, Y., Wang, Z.M., Tan, W., Wang, X., Li, Y., Bai, B., Li, Y., Zhang, S.F., Yan, H.L., Chen, Z.L. and Liu, C.M. (2018). Partial loss of psychiatric risk gene Mir137 in mice causes repetitive behavior and impairs sociability and learning via increased Pde10a. *Nature neuroscience*, 21(12):1689-1703.
53. Chin, E.W., Lim, W.M., Ma, D., Rosales, F.J. and Goh, E.L. (2018). Choline Rescues Behavioural Deficits in a Mouse Model of Rett Syndrome by Modulating Neuronal Plasticity. *Molecular neurobiology*, PubMed PMID: 30220058.
54. Choi, J.Y., Mohibbullah, M., Park, I.S., Moon, I.S. and Hong, Y.K. (2018). An ethanol extract from the phaeophyte *Undaria pinnatifida* improves learning and memory impairment and dendritic spine morphology in hippocampal neurons. *Journal of Applied Phycology*, 30(1):129-136.

55. Choi, S.H., Bylykbashi, E., Chatila, Z.K., Lee, S.W., Pulli, B., Clemenson, G.D., Kim, E., Rompala, A., Oram, M.K., Asselin, C. and Aronson, J., (2018). Combined adult neurogenesis and BDNF mimic exercise effects on cognition in an Alzheimer's mouse model. *Science*, 361(6406), p.eaan8821.
56. Cicvaric, A., Yang, J., Bulat, T., Zambon, A., Dominguez-Rodriguez, M., Kühn, R., Sadowicz, M.G., Siwert, A., Egea, J., Pollak, D.D. and Moeslinger, T. (2018). Enhanced synaptic plasticity and spatial memory in female but not male FLRT2-haplodeficient mice. *Scientific reports*, 8(1):3703.
57. Cohen, S.M., Suutari, B., He, X., Wang, Y., Sanchez, S., Tirko, N.N., Mandelberg, N.J., Mullins, C., Zhou, G., Wang, S. and Kats, I. (2018). Calmodulin shuttling mediates cytonuclear signaling to trigger experience-dependent transcription and memory. *Nature Communications*, 9(1):2451.
58. Dai, L.J., Huang, Y.L., Sun, W.W., Ye, B.F. and Chen, S.Q., (2018). Increased glutamic acid decarboxylase expression in the hippocampus impairs learning and memory in Fmr1 knockout mice. *International Journal of Clinical and Experimental Medicine*, 11(4), pp.3121-3134.
59. Davenport, E.C., Szulc, B., Drew, J., Taylor, J., Morgan, T., Higgs, N.F., Lopez-Domenech, G. and Kittler, J.T. (2018). Correct CYFIP1 dosage is essential for synaptic inhibition and the excitatory/inhibitory balance. *bioRxiv*, p.303446.
60. de la Parra, J., Cuartero, M.I., Pérez-Ruiz, A., García-Culebras, A., Martín, R., Sánchez-Prieto, J., García-Segura, J.M., Lizasoain, I. and Moro, M.A. (2018). AhR Deletion Promotes Aberrant Morphogenesis and Synaptic Activity of Adult-Generated Granule Neurons and Impairs Hippocampus-Dependent Memory. *eNeuro*, PubMed PMID: 30225360.
61. De Munter, S., Bamps, D., Malheiro, A.R., Kumar Baboota, R., Brites, P. and Baes, M. (2018). Autonomous Purkinje cell axonal dystrophy causes ataxia in peroxisomal multifunctional protein-2 deficiency. *Brain Pathology*, 28(5):631-643.
62. Dong, E., Guidotti, A., Zhang, H. and Pandey, S.C., (2018). Prenatal stress leads to chromatin and synaptic remodeling and excessive alcohol intake comorbid with anxiety-like behaviors in adult offspring. *Neuropharmacology*, 140, pp.76-85.
63. Dong, H., Keegan, J.M., Hong, E., Gallardo, C., Montalvo-Ortiz, J., Wang, B., Rice, K.C. and Csernansky, J. (2018). Corticotrophin releasing factor receptor 1 antagonists prevent chronic stress-induced behavioral changes and synapse loss in aged rats. *Psychoneuroendocrinology*, 90:92-101.
64. Dong, H., Zhu, M., Meng, L., Ding, Y., Yang, D., Zhang, S., Qiang, W., Fisher, D.W. and Xu, E.Y. (2018). Pumilio2 regulates synaptic plasticity via translational repression of synaptic receptors in mice. *Oncotarget*, 9(63):32134-32148.
65. Doria, J.G., de Souza, J.M., Silva, F.R., Olmo, I.G., Carvalho, T.G., Alves-Silva, J., Ferreira-Vieira, T.H., Santos, J.T., Xavier, C.Q., Silva, N.C. and Maciel, E.M. (2018). The mGluR5 positive allosteric modulator VU0409551 improves synaptic plasticity and memory of a mouse model of Huntington's disease. *Journal of neurochemistry*, 147(2):222-239.
66. Evanson, N.K., Guilhaume-Correa, F., Herman, J.P. and Goodman, M.D., (2018). Optic tract injury after closed head traumatic brain injury in mice: A model of indirect traumatic optic neuropathy. *PLoS one*, 13(5), p.e0197346.
67. Fan, C., Song, Q., Wang, P., Li, Y., Yang, M. and Yu, S.Y. (2018). Neuroprotective effects of ginsenoside-Rg1 against depression-like behaviors via suppressing glial activation, synaptic deficits and neuronal apoptosis in rats. *Frontiers in Immunology*, 9:2889.
68. Fan, C., Song, Q., Wang, P., Li, Y., Yang, M., Liu, B. and Yu, S.Y. (2018). Curcumin protects against chronic stress-induced dysregulation of neuroplasticity and depression-like behaviors via suppressing IL-1 β pathway in rats. *Neuroscience*, 392:92-106.
69. Fan, C., Zhu, X., Song, Q., Wang, P., Liu, Z. and Yu, S.Y. (2018). MiR-134 modulates chronic stress-induced structural plasticity and depression-like behaviors via downregulation of Limk1/cofilin signaling in rats. *Neuropharmacology*, 131:364-376.
70. Fogaça, M.V., Campos, A.C., Coelho, L.D., Duman, R.S. and Guimarães, F.S. (2018). The anxiolytic effects of cannabidiol in chronically stressed mice are mediated by the endocannabinoid system: role of neurogenesis and dendritic remodeling. *Neuropharmacology*, 135:22-33.
71. Frederick, N.M., Shah, P.V., Didonna, A., Langley, M.R., Kanthasamy, A.G. and Opal, P. (2018). Loss of the dystonia gene Thap1 leads to transcriptional deficits that converge on common pathogenic pathways in dystonic syndromes. *Human Molecular Genetics*, PubMed PMID: 30590536.
72. Fuchs, C., Fustini, N., Trazzi, S., Gennaccaro, L., Rimondini, R. and Ciani, E. (2018). Treatment with the GSK 3- β inhibitor Tideglusib improves hippocampal development and memory performance in juvenile, but not adult, Cdkl5 knockout mice. *European Journal of Neuroscience*, 47(9):1054-1066.
73. Fuchs, C., Gennaccaro, L., Trazzi, S., Bastianini, S., Bettini, S., Martire, V.L., Ren, E., Medici, G., Zoccoli, G., Rimondini, R. and Ciani, E. (2018). Heterozygous CDKL5 Knockout Female Mice Are a Valuable Animal Model for CDKL5 Disorder. *Neural Plasticity*, PubMed PMID: 29977282.
74. Garrett, L., Ung, M.C., Heermann, T., Niedermeier, K.M. and Hölter, S., (2018). Analysis of Neuropsychiatric Disease-Related Functional Neuroanatomical Markers in Mice. *Current Protocols in Mouse Biology*, 8(1), pp.79-128.
75. Goeke, C.M., Roberts, M.L., Hashimoto, J.G., Finn, D.A. and Guizzetti, M. (2018). Neonatal Ethanol and Choline Treatments Alter the Morphology of Developing Rat Hippocampal Pyramidal Neurons in Opposite

- Directions. *Neuroscience*, 374:13-24.
76. Gray, N.E., Zweig, J.A., Caruso, M., Martin, M.D., Zhu, J.Y., Quinn, J.F. and Soumyanath, A. (2018). Centella asiatica increases hippocampal synaptic density and improves memory and executive function in aged mice. *Brain and Behavior*, 8(7):e01024.
 77. Griffiths, K.K., Wang, A., Wang, L., Tracey, M., Kleiner, G., Quinzii, C.M., Perez-Zoghbi, J.F., Licznarski, P., Jonas, E.A. and Levy, R.J., (2018). Inefficient Thermogenic Mitochondrial Respiration Due to Futile Proton Leak in a Mouse Model of Fragile X Syndrome. Available at SSRN 3287782.
 78. Gross, C., Banerjee, A., Tiwari, D., Longo, F., White, A.R., Allen, A.G., Schroeder-Carter, L.M., Krzeski, J.C., Elsayed, N.A., Puckett, R. and Klann, E. (2018). Isoform-selective phosphoinositide 3-kinase inhibition ameliorates a broad range of fragile X syndrome-associated deficits in a mouse model. *Neuropsychopharmacology*, PubMed PMID: 30061744.
 79. Gstrein, T., Edwards, A., Přistoupilová, A., Leca, I., Breuss, M., Pilat-Carotta, S., Hansen, A.H., Tripathy, R., Traunbauer, A.K., Hochstoeger, T. and Rosoklija, G. (2018). Mutations in Vps15 perturb neuronal migration in mice and are associated with neurodevelopmental disease in humans. *Nature neuroscience*, 21(2):207.
 80. Guo, Y., Benson, C., Hill, M., Henry, S., Efraim, P., Waxman, S.G., Dib-Hajj, S. and Tan, A.M. (2018). Therapeutic potential of Pak1 inhibition for pain associated with cutaneous burn injury. *Molecular pain*, PubMed PMID: 29956587
 81. Hansen, S., Jørgensen, J., Nyegaard, J., Lykkesfeldt, J. and Tveden-Nyborg, P. (2018). Early life vitamin C deficiency does not alter morphology of hippocampal CA1 pyramidal neurons or markers of synaptic plasticity in a Guinea pig model. *Nutrients*, 10(6):749.
 82. Hayase, S., Wang, H., Ohgushi, E., Kobayashi, M., Mori, C., Horita, H., Mineta, K., Liu, W.C. and Wada, K. (2018). Vocal practice regulates singing activity-dependent genes underlying age-independent vocal learning in songbirds. *PLoS biology*, 16(9):e2006537.
 83. Herring, A. and Keyvani, K., Universitat Duisburg-Essen (2018). Agents inhibiting kallikrein-8 for use in the prevention or treatment of alzheimer's disease. *U.S. Patent Application* 15/761,725.
 84. Højland, A., Richner, M., Mølgaard, S., Dieu, R.S., Eskelund, A., Nykjer, A., Nyegaard, J.R., Lykkesfeldt, J., Glerup, S. and Nielsen, M.S. (2018). Biochemical and cognitive effects of docosahexaenoic acid differ in a developmental and SorLA dependent manner. *Behavioural brain research*, 348:90-100.
 85. Hongsheng, M., Li, R., Han, C., Lu, X. and Zhang, H. (2018). Minocycline promotes post-hemorrhagic neurogenesis via M2 microglia polarization via up-regulation of the TrkB/BDNF pathway in rats. *Journal of neurophysiology*, 120(3):1307-1317.
 86. Hosseini, S., Wilk, E., Michaelsen-Preusse, K., Gerhauser, I., Baumgärtner, W., Geffers, R., Schughart, K. and Korte, M. (2018). Long-term neuroinflammation induced by influenza A virus infection and the impact on hippocampal neuron morphology and function. *Journal of Neuroscience*, 38(12):3060-3080.
 87. Hottman, D., Cheng, S., Gram, A., LeBlanc, K., Yuan, L.L. and Li, L. (2018). Systemic or forebrain neuron-specific deficiency of geranylgeranyltransferase-1 impairs synaptic plasticity and reduces dendritic spine density. *Neuroscience*, 373:207-217.
 88. Huang, L., Shum, E.Y., Jones, S.H., Lou, C.H., Dumdie, J., Kim, H., Roberts, A.J., Jolly, L.A., Espinoza, J.L., Skarbrevik, D.M. and Phan, M.H., (2018). A Upf3b-mutant mouse model with behavioral and neurogenesis defects. *Molecular psychiatry*, 23(8), p.1773.
 89. Huang, S.H., Wu, W.R., Lee, L.M., Huang, P.R. and Chen, J.C. (2018). mTOR signaling in the nucleus accumbens mediates behavioral sensitization to methamphetamine. *Progress in Neuro-Psychopharmacology and Biological Psychiatry*, 86:331-339.
 90. Ignatenko, O., Chilov, D., Paetau, I., Miguel, E., Jackson, C.B., Capin, G., Paetau, A., Terzioglu, M., Euro, L. and Suomalainen, A. (2018). Loss of mtDNA activates astrocytes and leads to spongiform encephalopathy. *Nature communications*, 9(1):70.
 91. Jeanneteau, F., Barrère, C., Vos, M., De Vries, C.J., Rouillard, C., Levesque, D., Dromard, Y., Moisan, M.P., Duric, V., Franklin, T.C. and Duman, R.S. (2018). The stress-induced transcription factor NR4A1 adjusts mitochondrial function and synapse number in prefrontal cortex. *Journal of Neuroscience*, 38(6):1335-1350.
 92. Ji, M., Yuan, H., Yuan, S., Xia, J. and Yang, J. (2018). The p75 neurotrophin receptor might mediate sepsis-induced synaptic and cognitive impairments. *Behavioural brain research*, 347:339-349.
 93. Jie, J., Xu, X., Xia, J., Tu, Z., Guo, Y., Li, C., Zhang, X., Wang, H., Song, W. and Xie, P. (2018). Memory Impairment Induced by Borna Disease Virus 1 Infection is Associated with Reduced H3K9 Acetylation. *Cellular Physiology and Biochemistry*, 49(1):381-394.
 94. Joseph, A.T., Bhardwaj, S.K. and Srivastava, L.K. (2018). Role of prefrontal cortex anti-and pro-inflammatory cytokines in the development of abnormal behaviors induced by disconnection of the ventral hippocampus in neonate rats. *Frontiers in Behavioral Neuroscience*, 12:244.
 95. Jun, M., Xiaolong, Q., Chaojuan, Y., Ruiyuan, P., Shukun, W., Junbing, W., Li, H., Hong, C., Jinbo, C., Rong, W. and Yajin, L., (2018). Calhm2 governs astrocytic ATP releasing in the development of depression-like behaviors. *Molecular psychiatry*, 23:883-891
 96. Kang, S., Lee, S., Kim, J., Kim, J.C., Kim, S.H., Son, Y., Shin, T., Youn, B., Kim, J.S., Wang, H. and Yang, M. (2018). Chronic Treatment with Combined Chemotherapeutic Agents Affects Hippocampal Micromorphometry and Function in Mice, Independently of Neuroinflammation. *Experimental neurobiology*, 27(5):419-436.

97. Kasza, Á., (2018). Rodent models of Alzheimer's disease: Intracerebroventricular administration of β amyloid 1-42 oligomers into wild type rats and testing LA1011 drug candidate on app/ps1 transgenic mice (*Doctoral dissertation*, University of Szeged, Hungary).
98. Keil, K.P., Sethi, S., Wilson, M.D., Silverman, J.L., Pessah, I.N. and Lein, P.J. (2018). Genetic mutations in Ca^{2+} signaling alter dendrite morphology and social approach in juvenile mice. *Genes, Brain and Behavior*, 12:e12526.
99. Keyvani, K., Münster, Y., Kurapati, N.K., Rubach, S., Schönborn, A., Kocakavuk, E., Karout, M., Hammesfahr, P., Wang, Y.C., Hermann, D.M. and Teuber-Hanselmann, S. (2018). Higher levels of kallikrein-8 in female brain may increase the risk for Alzheimer's disease. *Brain Pathology*, 28(6):947-964.
100. Khatri, N., Gilbert, J.P., Huo, Y., Sharafrazi, R., Nee, M., Qiao, H. and Man, H.Y. (2018). The autism protein Ube3A/E6AP remodels neuronal dendritic arborization via caspase-dependent microtubule destabilization. *Journal of Neuroscience*, 38(2):363-378.
101. Klahr, A.C., Fagan, K., Aziz, J.R., John, R. and Colbourne, F., (2018). Mild Contralesional Hypothermia Reduces Use of the Unimpaired Forelimb in a Skilled Reaching Task After Motor Cortex Injury in Rats. *Therapeutic hypothermia and temperature management*, 8(2), pp.90-98.
102. Klok, M.D., Bugiani, M., de Vries, S.I., Gerritsen, W., Breur, M., van der Sluis, S., Heine, V.M., Kole, M.H., Baron, W. and van der Knaap, M.S. (2018). Axonal abnormalities in vanishing white matter. *Annals of clinical and translational neurology*, 5(4):429-444.
103. Kumamoto, H., Yamaguchi, T., Konno, K., Izumi, T., Yoshida, T., Ohmura, Y., Watanabe, M. and Yoshioka, M. (2018). Repeated fluvoxamine treatment recovers early postnatal stress-induced hypersociability-like behavior in adult rats. *Journal of pharmacological sciences*, 136(1):1-8.
104. Lan, Y., Sullivan, P.M. and Hu, F. (2018). SMCR8 negatively regulates AKT and MTORC1 signaling to modulate lysosome biogenesis and tissue homeostasis. *bioRxiv*, p.485243.
105. Lee, F.Y., Wang, H.B., Hitchcock, O.N., Loh, D.H., Whittaker, D.S., Kim, Y.S., Aiken, A., Kokikian, C., Dell'Angelica, E.C., Colwell, C.S. and Ghiani, C.A. (2018). Sleep/Wake Disruption in a Mouse Model of BLOC-1 Deficiency. *Frontiers in neuroscience*, 12:759.
106. Leng, L., Zhuang, K., Liu, Z., Huang, C., Gao, Y., Chen, G., Lin, H., Hu, Y., Wu, D., Shi, M. and Xie, W. (2018). Menin deficiency leads to depressive-like behaviors in mice by modulating astrocyte-mediated neuroinflammation. *Neuron*, 100(3):551-563.
107. Lewis, M.H., Lindenmaier, Z., Boswell, K., Edington, G., King, M.A. and Muehlmann, A.M. (2018). Subthalamic nucleus pathology contributes to repetitive behavior expression and is reversed by environmental enrichment. *Genes, Brain and Behavior*, 17(8):e12468.
108. Li, J., Zhang, S., Li, C., Li, M. and Ma, L. (2018). Sitagliptin rescues memory deficits in Parkinsonian rats via upregulating BDNF to prevent neuron and dendritic spine loss. *Neurological research*, 40(9):736-743.
109. Liu, L., Xuan, C., Shen, P., He, T., Chang, Y., Shi, L., Tao, S., Yu, Z., Brown, R.E. and Wang, J. (2018). Hippocampal Mechanisms Underlying Impairment in Spatial Learning Long After Establishment of Noise-Induced Hearing Loss in CBA Mice. *Frontiers in systems neuroscience*, 12:35.
110. Lu, L., Ben, X., Xiao, L., Peng, M. and Zhang, Y. (2018). AMP-activated protein kinase activation in mediating phenylalanine-induced neurotoxicity in experimental models of phenylketonuria. *Journal of Inherited Metabolic Disease: Official Journal of the Society for the Study of Inborn Errors of Metabolism*, 41(4):679-687.
111. Mahmood, U., Ahn, S., Yang, E.J., Choi, M., Kim, H., Regan, P., Cho, K. and Kim, H.S. (2018). Dendritic spine anomalies and PTEN alterations in a mouse model of VPA-induced autism spectrum disorder. *Pharmacological research*, 128:110-121.
112. Mancezak, M., Kandimalla, R., Yin, X. and Reddy, P.H., (2018). Hippocampal mutant APP and amyloid beta-induced cognitive decline, dendritic spine loss, defective autophagy, mitophagy and mitochondrial abnormalities in a mouse model of Alzheimer's disease. *Human molecular genetics*, 27(8), pp.1332-1342.
113. Martyn, A.C., Toth, K., Schmalzigaug, R., Hedrick, N.G., Rodriguez, R.M., Yasuda, R., Wetsel, W.C. and Premont, R.T. (2018). GIT1 regulates synaptic structural plasticity underlying learning. *PLoS one*, 13(3):e0194350.
114. Matagne, V., Wondolowski, J., Frerking, M., Shahidullah, M., Delamere, N.A., Sandau, U.S., Budden, S. and Ojeda, S.R., (2018). Correcting deregulated Fxyd1 expression rescues deficits in neuronal arborization and potassium homeostasis in MeCP2 deficient male mice. *Brain research*, 1697, pp.45-52.
115. Moore, A.R., Richards, S.E., Kenny, K., Royer, L., Chan, U., Flavahan, K., Van Hooser, S.D. and Paradis, S. (2018). Rem2 stabilizes intrinsic excitability and spontaneous firing in visual circuits. *Elife*, 7:e33092.
116. Morel, C., Sherrin, T., Kennedy, N.J., Forest, K.H., Barutcu, S.A., Robles, M., Carpenter-Hyland, E., Alfulajj, N., Standen, C.L., Nichols, R.A. and Benveniste, M. (2018). JIP1-Mediated JNK Activation Negatively Regulates Synaptic Plasticity and Spatial Memory. *Journal of Neuroscience*, 38(15):3708-3728.
117. Morud, J., Strandberg, J., Andrén, A., Ericson, M., Söderpalm, B. and Adermark, L. (2018). Progressive modulation of accumbal neurotransmission and anxiety-like behavior following protracted nicotine withdrawal. *Neuropharmacology*, 128:86-95.
118. Nagarajan, N., Jones, B.W., West, P.J., Marc, R.E. and Capecchi, M.R. (2018). Corticostriatal circuit defects in Hoxb8 mutant mice. *Molecular psychiatry*, 23(9):1.
119. Nagpal, L., Fu, C. and Snyder, S.H. (2018). Inositol Hexakisphosphate Kinase-2 in Cerebellar Granule Cells

- Regulates Purkinje Cells and Motor Coordination via Protein 4.1 N. *Journal of Neuroscience*, 38(34):7409-7419.
120. Nelissen, T.P., Bamford, R.A., Tochitani, S., Akkus, K., Kudzinskas, A., Yokoi, K., Okamoto, H., Yamamoto, Y., Burbach, J.P.H., Matsuzaki, H. and Oguro-Ando, A. (2018). CD38 is required for dendritic organization in visual cortex and hippocampus. *Neuroscience*, 372:114-125.
121. Nestor, J., Arinuma, Y., Huerta, T.S., Kowal, C., Nasiri, E., Kello, N., Fujieda, Y., Bialas, A., Hammond, T., Sriram, U. and Stevens, B. (2018). Lupus antibodies induce behavioral changes mediated by microglia and blocked by ACE inhibitors. *Journal of Experimental Medicine*, 215(10):2554-2566.
122. Nie, X., Kitaoka, S., Tanaka, K., Segi-Nishida, E., Imoto, Y., Ogawa, A., Nakano, F., Tomohiro, A., Nakayama, K., Taniguchi, M. and Mimori-Kiyosue, Y., (2018). The innate immune receptors TLR2/4 mediate repeated social defeat stress-induced social avoidance through prefrontal microglial activation. *Neuron*, 99(3), pp.464-479.
123. Nimmagadda, K. and Weiland, J.D. (2018). Retinotopic Responses in the Visual Cortex Elicited by Epiretinal Electrical Stimulation in Normal and Retinal Degenerate Rats. *Translational vision science & technology*, 7(5):33-33.
124. Novaes, L.S., dos Santos, N.B., Perfetto, J.G., Goosens, K.A. and Munhoz, C.D. (2018). Environmental enrichment prevents acute restraint stress-induced anxiety-related behavior but not changes in basolateral amygdala spine density. *Psychoneuroendocrinology*, 98:6-10.
125. Nowak, D., De Groef, L., Moons, L. and Mozrzymas, J.W. (2018). MMP-3 deficiency does not influence the length and number of CA1 dendrites of hippocampus of adult mice. *Acta neurobiologiae experimentalis*, 78(3):281-286.
126. Nuber, S., Rajsombath, M., Minakaki, G., Winkler, J., Müller, C.P., Ericsson, M., Caldarone, B., Dettmer, U. and Selkoe, D.J., (2018). Abrogating native α -synuclein tetramers in mice causes a L-DOPA-responsive motor syndrome closely resembling Parkinson's disease. *Neuron*, 100(1), pp.75-90.
127. O'Reilly, K.C., Levy, E.R., Patino, A.V., Perica, M.I. and Fenton, A.A. (2018). Sub-circuit alterations in dorsal hippocampus structure and function after global neurodevelopmental insult. *Brain Structure and Function*, 223(8):3543-3556.
128. Ogundeke, O.M. and Lee, C.C., (2018). CaMKII α expression in a mouse model of NMDAR hypofunction schizophrenia: Putative roles for IGF-1R and TLR4. *Brain research bulletin*, 137, pp.53-70.
129. Okuda, K., Takao, K., Watanabe, A., Miyakawa, T., Mizuguchi, M. and Tanaka, T. (2018). Comprehensive behavioral analysis of the Cdk5 knockout mice revealed significant enhancement in anxiety-and fear-related behaviors and impairment in both acquisition and long-term retention of spatial reference memory. *PLoS one*, 13(4):e0196587.
130. Olivares-Hernández, J.D., García-García, F., Camacho-Abrego, I., Flores, G. and Juárez-Aguilar, E. (2018). Intracerebroventricular administration of growth hormone induces morphological changes in pyramidal neurons of the hippocampus and prefrontal cortex in adult rats. *Synapse*, 72(7):e22030.
131. Ortiz, J.B., Anglin, J.M., Daas, E.J., Paode, P.R., Nishimura, K. and Conrad, C.D. (2018). BDNF and TrkB Mediate the Improvement from Chronic Stress-induced Spatial Memory Deficits and CA3 Dendritic Retraction. *Neuroscience*, 388:330-346.
132. Paramo, B., Wyatt, S. and Davies, A.M. (2018). An essential role for neuregulin-4 in the growth and elaboration of developing neocortical pyramidal dendrites. *Experimental neurology*, 302:85-92.
133. Park, M.H., Lee, J.Y., Park, K.H., Jung, I.K., Kim, K.T., Lee, Y.S., Ryu, H.H., Jeong, Y., Kang, M., Schwanger, M. and Gulbins, E., (2018). Vascular and neurogenic rejuvenation in aging mice by modulation of ASM. *Neuron*, 100(1), pp.167-182.
134. Parras, A., Anta, H., Santos-Galindo, M., Swarup, V., Elorza, A., Nieto-González, J.L., Picó, S., Hernández, I.H., Díaz-Hernández, J.I., Belloc, E. and Rodolose, A., (2018). Autism-like phenotype and risk gene mRNA deadenylation by CPEB4 mis-splicing. *Nature*, 560(7719), p.441.
135. Paternoster, V., Rajkumar, A.P., Nyengaard, J.R., Borglum, A.D., Grove, J. and Christensen, J.H. (2018). The importance of data structure in statistical analysis of dendritic spine morphology. *Journal of neuroscience methods*, 296:93-98.
136. Pearn, M.L., Schilling, J.M., Jian, M., Egawa, J., Wu, C., Mandyam, C.D., Fannon-Pavlich, M.J., Nguyen, U., Bertoglio, J., Kodama, M. and Mahata, S.K., (2018). Inhibition of RhoA reduces propofol-mediated growth cone collapse, axonal transport impairment, loss of synaptic connectivity, and behavioural deficits. *British journal of anaesthesia*, 120(4), pp.745-760.
137. Peng, B., Peng, S.H., Qu, R.M., Xu, L.H. and Jiang, Z.L. (2018). Nitrogen narcosis induced by repetitive hyperbaric nitrogen oxygen mixture exposure impairs long-term cognitive function in newborn mice. *PLoS one*, 13(4):e0196611.
138. Pérez, M.Á., Peñalosa-Sancho, V., Ahumada, J., Fuenzalida, M. and Dagnino-Subiabre, A. (2018). n-3 Polyunsaturated fatty acid supplementation restored impaired memory and GABAergic synaptic efficacy in the hippocampus of stressed rats. *Nutritional neuroscience*, 21(8):556-569.
139. Petit-Pedrol, M., Sell, J., Planagumà, J., Mannara, F., Radosevic, M., Haselmann, H., Ceanga, M., Sabater, L., Spatola, M., Soto, D. and Gasull, X., (2018). LGI1 antibodies alter Kv1.1 and AMPA receptors changing synaptic excitability, plasticity and memory. *Brain*, 141(11), pp.3144-3159.

140. Pratchayasakul, W., Thongnak, L.O., Chattipakorn, K., Lungaphin, A., Pongchaidecha, A., Satjaritanun, P., Jaiwongkam, T., Kerdphoo, S. and Chattipakorn, S.C. (2018). Atorvastatin and insulin equally mitigate brain pathology in diabetic rats. *Toxicology and applied pharmacology*, 342:79-85.
141. Qu, Y., Yang, C., Ren, Q., Ma, M., Dong, C. and Hashimoto, K. (2018). Regional differences in dendritic spine density confer resilience to chronic social defeat stress. *Acta neuropsychiatrica*, 30(2):117-122.
142. Quach, T., Auvergnon, N., Khanna, R., Belin, M.F., Kolattukudy, P., Honnorat, J. and Duchemin, A.M. (2018). Opposing Morphogenetic Defects on Dendrites and Mossy Fibers of Dentate Granular Neurons in CRMP3-Deficient Mice. *Brain sciences*, 8(11):196.
143. Rajkumar, A.P., Qvist, P., Larsen, S.H., Lazarus, R., Pallesen, J., Nava, N., Winther, G., Liebenberg, N., Paternoster, V., Fryland, T. and Palmfeldt, J. (2018). The neurobiology of BRD1 implicates sex-biased dysregulation of nuclear receptor signaling in mental disorders. *bioRxiv*, p.257170.
144. Ramos-Fernández, E., Tapia-Rojas, C., Ramirez, V.T. and Inestrosa, N.C. (2018). Wnt-7a Stimulates Dendritic Spine Morphogenesis and PSD-95 Expression Through Canonical Signaling. *Molecular neurobiology*, PubMed PMID: 29967987.
145. Ran, Y.H., Hu, X.X., Wang, Y.L., Zhao, N., Zhang, L.M., Liu, H.X. and Li, Y.F. (2018). YL-0919, a dual 5-HT_{1A} partial agonist and SSRI, produces antidepressant and anxiolytic-like effects in rats subjected to chronic unpredictable stress. *Acta Pharmacologica Sinica*, 39(1):12.
146. Raven, F., Meerlo, P., Van der Zee, E.A., Abel, T. and Havekes, R. (2018). A brief period of sleep deprivation causes spine loss in the dentate gyrus of mice. *Neurobiology of learning and memory*, PubMed PMID: 29588221.
147. Reeber, S.L., Arancillo, M. and Sillitoe, R.V. (2018). Bergmann glia are patterned into topographic molecular zones in the developing and adult mouse cerebellum. *The Cerebellum*, 17(4):392-403.
148. Reker, A.N., Oliveros, A., Sullivan, J.M., Nahar, L., Hinton, D.J., Kim, T., Bruner, R.C., Choi, D.S., Goeders, N.E. and Nam, H.W. (2018). Neurogranin in the nucleus accumbens regulates NMDA receptor tolerance and motivation for ethanol seeking. *Neuropharmacology*, 131:58-67.
149. Ren, Q.G., Gong, W.G., Zhou, H., Shu, H., Wang, Y.J. and Zhang, Z.J. (2018). Spatial Training Ameliorates Long-Term Alzheimer's Disease-Like Pathological Deficits by Reducing NLRP3 Inflammasomes in PR5 Mice. *Neurotherapeutics*, PubMed PMID: 3056048.
150. Rincel, M., Lépinay, A.L., Janthakhin, Y., Soudain, G., Yvon, S., Da Silva, S., Joffre, C., Aubert, A., Séré, A., Layé, S. and Theodorou, V. (2018). Maternal high-fat diet and early life stress differentially modulate spine density and dendritic morphology in the medial prefrontal cortex of juvenile and adult rats. *Brain Structure and Function*, 223(2):883-895.
151. Risher, W.C., Kim, N., Koh, S., Choi, J.E., Mitev, P., Spence, E.F., Pilaz, L.J., Wang, D., Feng, G., Silver, D.L. and Soderling, S.H. (2018). Thrombospondin receptor $\alpha 2\delta$ -1 promotes synaptogenesis and spinogenesis via postsynaptic Rac1. *J Cell Biology*, 217(10):3747-3765.
152. Romain, F.O.L., Braudeau, J., Cartier, N., Buchholz, C., Tobias, A. and Mueller, U., (2018). Methods and pharmaceutical composition for the treatment of alzheimer's disease. *U.S. Patent Application* 15/580,934.
153. Rousseaux, M.W., Tschumperlin, T., Lu, H.C., Lackey, E.P., Bondar, V.V., Wan, Y.W., Tan, Q., Adamski, C.J., Friedrich, J., Twaroski, K. and Chen, W. (2018). ATXN1-C1C Complex Is the Primary Driver of Cerebellar Pathology in Spinocerebellar Ataxia Type 1 through a Gain-of-Function Mechanism. *Neuron*, 97(6):1235-1243.
154. Saito, Y., Yuan, Y., Zucker-Scharff, I., Fak, J.J., Jereb, S., Tajima, Y., Licatalosi, D.D. and Darnell, R.B., (2018). Differential NOVA2-Mediated Splicing in Excitatory and Inhibitory Neurons Regulates Cortical Development and Cerebellar Function. *Neuron*. PMID: 30638744.
155. Saito, Y., Yuan, Y., Zucker-Scharff, I., Fak, J.J., Tajima, Y., Licatalosi, D.D. and Darnell, R.B. (2018). Neuron-specific cTag-CLIP reveals cell-specific diversity of functional RNA regulation in the brain. *bioRxiv*, p.244905.
156. Sales, A.J., Fogaca, M.V., Sartim, A.G., Pereira, V.S., Wegener, G., Guimarães, F.S. and Joca, S.R. (2018). Cannabidiol Induces Rapid and Sustained Antidepressant-Like Effects Through Increased BDNF Signaling and Synaptogenesis in the Prefrontal Cortex. *Molecular neurobiology*, PubMed PMID: 29869197.
157. Sa-nguanmoo, P., Tanajak, P., Kerdphoo, S., Jaiwongkam, T., Wang, X., Liang, G., Li, X., Jiang, C., Pratchayasakul, W., Chattipakorn, N. and Chattipakorn, S.C. (2018). FGF21 and DPP-4 inhibitor equally prevents cognitive decline in obese rats. *Biomedicine & Pharmacotherapy*, 97:1663-1672.
158. Scarante, F.F., (2018). Behavioral and neuroplastic effects of the combination of escitalopram and cannabinoids (*Doctoral dissertation*, University of São Paulo, Brazil).
159. Scheuer, T., Sharkovska, Y., Tarabykin, V., Marggraf, K., Brockmöller, V., Bührer, C., Endesfelder, S. and Schmitz, T. (2018). Neonatal hyperoxia perturbs neuronal development in the cerebellum. *Molecular neurobiology*, 55(5):3901-3915.
160. Schick, A. (2018). *Overexpression of transmembrane-agrin in the murine central nervous system (Doctoral dissertation*, Ludwig-Maximilians-Universität zu München).
161. Shen, Y., Zhao, H., Li, P., Peng, Y., Cui, P., Miao, F., Zhang, Y., Zhang, A. and Zhang, J. (2018). MHC Class I Molecules and PirB Shape Neuronal Morphology by Affecting the Dendritic Arborization of Cortical Neurons. *Neurochemical research*, PubMed PMID: 30406910.
162. Shin, B.C., Cepeda, C., Estrada-Sánchez, A.M., Levine, M.S., Hodaei, L., Dai, Y., Jung, J., Ganguly, A., Clark, P. and Devaskar, S.U. (2018). Neural Deletion of Glucose Transporter Isoform 3 Creates Distinct Postnatal and

- Adult Neurobehavioral Phenotypes. *Journal of Neuroscience*, 38(44):9579-9599.
163. Smith, L.A. and McMahon, L.L. (2018). Deficits in synaptic function occur at medial perforant path-dentate granule cell synapses prior to Schaffer collateral-CA1 pyramidal cell synapses in the novel TgF344-Alzheimer's Disease Rat Model. *Neurobiology of disease*, 110:166-179.
 164. Sobue, A., Ito, N., Nagai, T., Shan, W., Hada, K., Nakajima, A., Murakami, Y., Mouri, A., Yamamoto, Y., Nabeshima, T. and Saito, K. (2018). Astroglial major histocompatibility complex class I following immune activation leads to behavioral and neuropathological changes. *Glia*, 66(5):1034-1052.
 165. Sobue, A., Kushima, I., Nagai, T., Shan, W., Kohno, T., Aleksic, B., Aoyama, Y., Mori, D., Arioka, Y., Kawano, N. and Yamamoto, M. (2018). Genetic and animal model analyses reveal the pathogenic role of a novel deletion of RELN in schizophrenia. *Scientific reports*, 8(1):13046
 166. Soteros, B.M., Cong, Q., Palmer, C.R. and Sia, G.M. (2018). Sociability and synapse subtype-specific defects in mice lacking SRPX2, a language-associated gene. *PLoS one*, 13(6):e0199399.
 167. Stanelle-Bertram, S., Walendy-Gnirß, K., Speiseder, T., Thiele, S., Asante, I.A., Dreier, C., Kouassi, N.M., Preuß, A., Pilitz-Stolze, G., Müller, U. and Thanisch, S. (2018). Male offspring born to mildly ZIKV-infected mice are at risk of developing neurocognitive disorders in adulthood. *Nature microbiology*, 3(10):1161.
 168. Suh, J., Romano, D.M., Herrick, S.P., Hooli, B., Oram, M.K., Mullin, K., Wasco, W., Schmammann, J.D., Albers, M.W. and Tanzi, R.E. (2018). Loss of Ataxin-1 Elevates BACE1 Expression and Impairs Axonal Targeting in the Cerebrum. Available at SSRN.3155885.
 169. Sun, J., Liu, Y., Jia, Y., Hao, X., Ju Lin, W., Tran, J., Lynch, G., Baudry, M. and Bi, X. (2018). UBE3A-mediated p18/LAMTOR1 ubiquitination and degradation regulate mTORC1 activity and synaptic plasticity. *Elife*, 7:e37993.
 170. Sun, M., Ahmad, N., Zhang, R. and Graw, J. (2018). Crybb2 associates with Tmsb4X and is crucial for dendrite morphogenesis. *Biochemical and biophysical research communications*, 503(1):123-130.
 171. Sun, X.D., Chen, W.B., Sun, D., Huang, J., Li, Y.Q., Pan, J.X., Wang, Y.N., Zhao, K., Dong, Z.Q., Wang, H.S. and Xiong, L. (2018). Neogenin in amygdala for neuronal activity and information processing. *Journal of Neuroscience*, 38(44):9600-9613.
 172. Synowiec, S., Lu, J., Yu, L., Goussakov, I., Lieber, R. and Drobyshvsky, A. (2018). Spinal Hyper-Excitability and Altered Muscle Structure Contribute to Muscle Hypertonia in Newborns After Antenatal Hypoxia-Ischemia in a Rabbit Cerebral Palsy Model. *Frontiers in Neurology*, 9:1183.
 173. Tang, J., Yu, W., Chen, S., Gao, Z. and Xiao, B. (2018). Microglia Polarization and Endoplasmic Reticulum Stress in Chronic Social Defeat Stress Induced Depression Mouse. *Neurochemical research*, 43(5):985-994.
 174. Tebbenkamp, A.T., Varela, L., Choi, J., Paredes, M.I., Giani, A.M., Song, J.E., Sestan-Pesa, M., Franjic, D., Sousa, A.M., Liu, Z.W. and Li, M. (2018). The 7q11.23 Protein DNAJC30 Interacts with ATP Synthase and Links Mitochondria to Brain Development. *Cell*, 175(4), pp.1088-1104.
 175. Thomas, T.C., Ogle, S.B., Rumney, B.M., May, H.G., Adelson, P.D. and Lifshitz, J. (2018). Does time heal all wounds? Experimental diffuse traumatic brain injury results in persisting histopathology in the thalamus. *Behavioral brain research*, 340:137-146.
 176. Trazzi, S., De Franceschi, M., Fuchs, C., Bastianini, S., Viggiano, R., Lupori, L., Mazziotti, R., Medici, G., Lo Martire, V., Ren, E. and Rimondini, R. (2018). CDKL5 protein substitution therapy rescues neurological phenotypes of a mouse model of CDKL5 disorder. *Human molecular genetics*, 27(9), pp.1572-1592.
 177. Uddin, M., Unda, B.K., Kwan, V., Holzapfel, N.T., White, S.H., Chalil, L., Woodbury-Smith, M., Ho, K.S., Harward, E., Murtaza, N. and Dave, B. (2018). OTUD7A regulates neurodevelopmental phenotypes in the 15q13.3 microdeletion syndrome. *The American Journal of Human Genetics*, 102(2):278-295.
 178. Ueda, Y., Bando, Y., Misumi, S., Ogawa, S., Ishida, A., Jung, C.G., Shimizu, T. and Hida, H. (2018). Alterations of Both Dendrite Morphology and Weaker Electrical Responsiveness in the Cortex of Hip Area Occur Before Rearrangement of the Motor Map in Neonatal White Matter Injury Model. *Frontiers in neurology*, 9. PMID: 29971036
 179. Ullman, J.C., Yang, J., Sullivan, M., Bendor, J., Levy, J., Pham, E., Silm, K., Seifkar, H., Sohal, V.S., Nicoll, R.A. and Edwards, R.H. (2018). A mouse model of autism implicates endosome pH in the regulation of pre-synaptic calcium entry. *Nature communications*, 9(1), p.330.
 180. van Campen, J.S., Hessel, E.V., Bohmbach, K., Rizzi, G., Lucassen, P.J., Lakshmi Turimella, S., Umeoka, E.H., Meerhoff, G.F., Braun, K.P., de Graan, P.N. and Joëls, M. (2018). stress and corticosteroids aggravate Morphological changes in the Dentate gyrus after early-life experimental Febrile seizures in Mice. *Frontiers in endocrinology*, 9:3.
 181. van Deijk, A.L.F. (2018). Astrocyte lipid metabolism in neuronal plasticity towards dietary modulation of synaptic function. (Doctoral dissertation, University in Amsterdam, Netherlands)
 182. Velazquez, R., Ferreira, E., Tran, A., Turner, E.C., Belfiore, R., Branca, C. and Oddo, S. (2018). Acute tau knockdown in the hippocampus of adult mice causes learning and memory deficits. *Aging cell*, 17(4):e12775.
 183. Vellema, M., Rocha, M., Bascones, S., Zsebok, S., Dreier, J., Leitner, S., Van der Linden, A., Brewer, J. and Gahr, M. (2018). Vocal motor experiences consolidate the vocal motor circuitry and accelerate future vocal skill development. *bioRxiv*, p.440388.
 184. Wahl, D., Solon-Biet, S.M., Wang, Q.P., Wali, J.A., Pulpitel, T., Clark, X., Raubenheimer, D., Senior, A.M., Sinclair, D.A., Cooney, G.J. and de Cabo, R. (2018). Comparing the Effects of Low-Protein and High-

- Carbohydrate Diets and Caloric Restriction on Brain Aging in Mice. *Cell reports*, 25(8):2234-2243.
185. Wang, H., Chen, F., Du, Y.F., Long, Y., Reed, M.N., Hu, M., Suppiramaniam, V., Hong, H. and Tang, S.S. (2018). Targeted inhibition of RAGE reduces amyloid- β influx across the blood-brain barrier and improves cognitive deficits in db/db mice. *Neuropharmacology*, 131:143-153.
186. Wang, J., Kou, X.L., Chen, C., Wang, M., Qi, C., Wang, J., You, W.Y., Hu, G., Chen, J. and Gao, J. (2018). Hippocampal Wdr1 Deficit Impairs Learning and Memory by Perturbing F-actin Depolymerization in Mice. *Cerebral Cortex*, PubMed PMID: 30590446.
187. Wang, R., Tan, J., Guo, J., Zheng, Y., Han, Q., So, K.F., Yu, J. and Zhang, L., (2018). Aberrant development and Wang, R., Tan, J., Guo, J., Zheng, Y., Han, Q., So, K.F., Yu, J. and Zhang, L., (2018). Aberrant development and synaptic transmission of cerebellar cortex in a VPA induced mouse autism model. *Frontiers in cellular neuroscience*, 12:500.
188. Wang, X., Li, R., Zacharek, A., Landschoot-Ward, J., Wang, F., Wu, K.H., Chopp, M., Chen, J. and Cui, X. (2018). Administration of Downstream ApoE Attenuates the Adverse Effect of Brain ABCA1 Deficiency on Stroke. *International journal of molecular sciences*, 19(11):3368.
189. Wang, X.Y., Zhou, H.R., Wang, S., Liu, C.Y., Qin, G.C., Fu, Q.Q., Zhou, J.Y. and Chen, L.X. (2018). NR2B-Tyr phosphorylation regulates synaptic plasticity in central sensitization in a chronic migraine rat model. *The journal of headache and pain*, 19(1):102.
190. Wang, Y., Zeng, C., Li, J., Zhou, Z., Ju, X., Xia, S., Li, Y., Liu, A., Teng, H., Zhang, K. and Shi, L. (2018). PAK2 Haploinsufficiency Results in Synaptic Cytoskeleton Impairment and Autism-Related Behavior. *Cell reports*, 24(8):2029-2041.
191. Widman, A.J., Stewart, A., Erb, E.M., Gardner, E. and McMahon, L.L. (2018). Intravascular ketamine increases theta-burst but not high frequency tetanus induced LTP at CA3-CA1 synapses within three hours and devoid of an increase in spine density. *Frontiers in synaptic neuroscience*, 10:8.
192. Wu, B., Li, C. and Lei, H. (2018). SAPAP4 Deletion Causes Synaptic Dysfunction in the nucleus accumbens. *Biochemical and biophysical research communications*, 505(4):1223-1227.
193. Wu, H. F., Chen, P. S., Hsu, Y. T., Lee, C. W., Wang, T. F., Chen, Y. J., & Lin, H. C. (2018). D-cycloserine ameliorates autism-like deficits by removing GluA2-containing AMPA receptors in a valproic acid-induced rat model. *Molecular Neurobiology*, 55:4811-24.
194. Wu, X., Lv, Y.G., Du, Y.F., Chen, F., Reed, M.N., Hu, M., Suppiramaniam, V., Tang, S.S. and Hong, H. (2018). Neuroprotective effects of INT-777 against A β 1-42-induced cognitive impairment, neuroinflammation, apoptosis, and synaptic dysfunction in mice. *Brain, Behavior, and Immunity*, 73:533-545.
195. Wu, Z., Li, X., Zhang, Y., Tong, D., Wang, L. and Zhao, P. (2018). Effects of sevoflurane exposure during mid-pregnancy on learning and memory in offspring rats: beneficial effects of maternal exercise. *Frontiers in cellular neuroscience*, 12:122.
196. Xiang, Q., Li, X.H., Fang, X.X., Jia, J., Ren, J., Dong, Y.C. and Ou-Yang, C. (2018). Alterations of synaptic plasticity in aged rats: Evidence of functional and morphological studies. *Technology and Health Care*, 26(2):271-278.
197. Xing, B., Han, G., Wang, M.J., Snyder, M.A. and Gao, W.J. (2018). Juvenile treatment with mGluR2/3 agonist prevents schizophrenia-like phenotypes in adult by acting through GSK3 β . *Neuropharmacology*, 137:359-371.
198. Xing, F.Z., Zhao, Y.G., Zhang, Y.Y., He, L., Zhao, J.K., Liu, M.Y., Liu, Y. and Zhang, J.Q. (2018). Nuclear and membrane estrogen receptor antagonists induce similar mTORC 2 activation-reversible changes in synaptic protein expression and actin polymerization in the mouse hippocampus. *CNS neuroscience & therapeutics*, 24(6):495-507.
199. Xu, Y., Cui, S.Y., Quan Ma, J.S., Yu, Y., Li, J.X., Zheng, L., Zhang, Y., Si, J.M. and Yu, Y.C. (2018). trans-Resveratrol ameliorates stress-induced irritable bowel syndrome-like behaviors by regulation of brain-gut axis. *Frontiers in pharmacology*, 9:63.
200. Yan, J., Porch, M.W., Bennett, M.V. and Zúkin, R.S., (2018). Activation of autophagy rescues synaptic and cognitive deficits in fragile X mice. *Proceedings of the National Academy of Sciences*, 115(41), pp.E9707-E9716.
201. Yang Eun Jung (2018). *A study on the effects of phloroglucinol on the cognitive impairments and neuropathological changes in Alzheimer's disease model* (Doctoral dissertation, Seoul National University Graduate School).
202. Yang, C., Kobayashi, S., Nakao, K., Dong, C., Han, M., Qu, Y., Ren, Q., Zhang, J.C., Ma, M., Toki, H. and Yamaguchi, J.I. (2018). AMPA receptor activation-independent antidepressant actions of ketamine metabolite (S)-norketamine. *Biological Psychiatry*, 84(8):591-600.
203. Yang, E.J., Mahmood, U., Kim, H., Choi, M., Choi, Y., Lee, J.P., Cho, J.Y., Hyun, J.W., Kim, Y.S., Chang, M.J. and Kim, H.S. (2018). Phloroglucinol ameliorates cognitive impairments by reducing the amyloid β peptide burden and pro-inflammatory cytokines in the hippocampus of 5XFAD mice. *Free Radical Biology and Medicine*, 126:221-234.
204. Yang, R., Walder-Christensen, K.K., Kim, N., Wu, D., Lorenzo, D., Badea, A., Gu, S., Streff, H., Soler-Alfonso, C., Meng, L. and Wetsel, W.C., (2018). A giant ankyrin-B mechanism for neuro-diversity/divergence through stochastic ectopic axon projections. *bioRxiv*, p.479741.
205. Yang, X., Cao, Z., Zhang, J., Shao, B., Song, M., Han, Y. and Li, Y. (2018). Dendritic spine loss caused by

- AIC13 is associated with inhibition of the Rac 1/cofilin signaling pathway. *Environmental Pollution*, 243:1689-1695.
206. Yin, J., Chen, W., Chao, E.S., Soriano, S., Wang, L., Wang, W., Cummock, S.E., Tao, H., Pang, K., Liu, Z. and Pereira, F.A. (2018). Otud7a knockout mice recapitulate many neurological features of 15q13.3 microdeletion syndrome. *The American Journal of Human Genetics*, 102(2):296-308.
207. Yoon, G., Shah, S.A., Ali, T. and Kim, M.O. (2018). The Adiponectin Homolog Osmotin Enhances Neurite Outgrowth and Synaptic Complexity via AdipoR1/NgR1 Signaling in Alzheimer's Disease. *Molecular neurobiology*, 55(8):6673-6686.
208. Yu, H., Zhong, J., Niu, B., Zhong, Q., Xiao, J., Xie, J., Lin, M., Zhou, Z., Xu, J. and Wang, H. (2018). Inhibition of Phosphodiesterase 4 by FCPR03 Alleviates Chronic Unpredictable Mild Stress-induced Depressive-like Behaviors and Prevents Dendritic Spine Loss in Mice Hippocampi. *International Journal of Neuropsychopharmacology*, PubMed PMID: 30407503.
209. Yu, Q., Liu, H., Sang, S., Chen, L., Zhao, Y., Wang, Y. and Zhong, C. (2018). Thiamine deficiency contributes to synapse and neural circuit defects. *Biological research*, 51(1):35.
210. Yu, Z., Wu, Y.J., Wang, Y.Z., Liu, D.S., Song, X.L., Jiang, Q., Li, Y., Zhang, S., Xu, N.J., Zhu, M.X. and Li, W.G., (2018). The acid-sensing ion channel ASIC1a mediates striatal synapse remodeling and procedural motor learning. *Sci. Signal.*, 11(542), p.eaar4481.
211. Zhang, J.H., Yu, L.J., Yang, H., Hui, Z., Jiang, S., Chen, L., Zhao, Y., Wang, S.L., Liu, Y. and Xu, Y. (2018). Huatuo Zaizao pill ameliorates cognitive impairment of APP/PS1 transgenic mice by improving synaptic plasticity and reducing A β deposition. *BMC complementary and alternative medicine*, 18(1):167.
212. Zhang, T., Jiang, X., Xu, M., Wang, H., Sang, X., Qin, M., Bao, P., Wang, R., Zhang, C., Lu, H. and Li, Y. (2018). Sleep and circadian abnormalities precede cognitive deficits in R521C FUS knockin rats. *Neurobiology of aging*, 72:159-170.
213. Zhang, W.J., Wang, H.H., Lv, Y.D., Liu, C.C., Sun, W.Y. and Tian, L.J. (2018). Downregulation of Egr-1 Expression Level via GluN2B Underlies the Antidepressant Effects of Ketamine in a Chronic Unpredictable Stress Animal Model of Depression. *Neuroscience*, 372:38-45.
214. Zhang, X. and Guizzetti, M. (2018). In Vitro and In Vitro Models for Alcohol Toxicity to the Extracellular Matrix of Astrocytes. In: *Methods in Pharmacology and Toxicology*. Humana Press.
215. Zhang, Y., Zhu, Z., Liang, H.Y., Zhang, L., Zhou, Q.G., Ni, H.Y., Luo, C.X. and Zhu, D.Y. (2018). nNOS-CAPON interaction mediates amyloid- β -induced neurotoxicity, especially in the early stages. *Aging cell*, 17(3):e12754.
216. Zhao, D., Meng, J., Zhao, Y., Huo, Y., Liu, Y., Zheng, N., Zhang, M., Gao, Y., Chen, Z., Sun, H. and Wang, X. (2018). RPS23RG1 Is Required for Synaptic Integrity and Rescues Alzheimer's Disease-Associated Cognitive Deficits. *Biological psychiatry*, PubMed PMID: 30292394.
217. Zhao, J., Bian, C., Liu, M., Zhao, Y., Sun, T., Xing, F. and Zhang, J. (2018). Orchiectomy and letrozole differentially regulate synaptic plasticity and spatial memory in a manner that is mediated by SRC-1 in the hippocampus of male mice. *The Journal of steroid biochemistry and molecular biology*, 178:354-368.
218. Zhao, J., Huynh, J., Hylin, M.J., O'Malley, J.J., Perez, A., Moore, A.N. and Dash, P.K. (2018). Mild traumatic brain injury reduces spine density of projection neurons in the medial prefrontal cortex and impairs extinction of contextual fear memory. *Journal of neurotrauma*, 35(1):149-156.
219. Zhao, Y., Wei, X., Song, J., Zhang, M., Huang, T. and Qin, J. (2018). Peroxisome Proliferator-Activated Receptor γ Agonist Rosiglitazone Protects Blood-Brain Barrier Integrity Following Diffuse Axonal Injury by Decreasing the Levels of Inflammatory Mediators Through a Caveolin-1-Dependent Pathway. *Inflammation*, PubMed PMID: 30488141.
220. Zhi, W.J., Peng, R.Y., Li, H.J., Zou, Y., Yao, B.W., Wang, C.Z., Liu, Z.H., Gao, X.H., Xu, X.P., Dong, J. and Zhao, L., (2018). Microwave radiation leading to shrinkage of dendritic spines in hippocampal neurons mediated by SNK-SPAR pathway. *Brain research*, 1679, pp.134-143.
221. Zhou, M., Gresack, J., Cheng, J., Uryu, K., Brichta, L., Greengard, P. and Flajolet, M. (2018). CK1 δ over-expressing mice display ADHD-like behaviors, frontostriatal neuronal abnormalities and altered expressions of ADHD-candidate genes. *Molecular Psychiatry*, 23: p.1.
222. Zhou, X., Xiao, W., Su, Z., Cheng, J., Zheng, C., Zhang, Z., Wang, Y., Wang, L., Xu, B., Li, S. and Yang, X., (2018). Hippocampal proteomic alteration in triple transgenic mouse model of Alzheimer's disease and implication of PINK 1 regulation in donepezil treatment. *Journal of proteome research*. PMID: 30484658
223. Zhu, B., Zhao, L., Luo, D., Xu, D., Tan, T., Dong, Z., Tang, Y., Min, Z., Deng, X., Sun, F. and Yan, Z. (2018). Furin promotes dendritic morphogenesis and learning and memory in transgenic mice. *Cellular and Molecular Life Sciences*, 75(13):2473-2488.
224. Zhu, C., Xu, Q., Mao, Z. and Lin, N., (2018). The Chinese Medicine Wu-Tou Decoction Relieves Neuropathic Pain by Inhibiting Hippocampal Microglia Activation. *Scientific reports*, 8(1), p.12292.
225. Zhu, H., Zhang, W., Zhao, Y., Shu, X., Wang, W., Wang, D., Yang, Y., He, Z., Wang, X. and Ying, Y. (2018). GSK3 β -mediated tau hyperphosphorylation triggers diabetic retinal neurodegeneration by disrupting synaptic and mitochondrial functions. *Molecular neurodegeneration*, 13(1):62.
226. Zhu, L.J., Ni, H.Y., Chen, R., Chang, L., Shi, H.J., Qiu, D., Zhang, Z., Wu, D.L., Jiang, Z.C., Xin, H.L. and Zhou, Q.G. (2018). Hippocampal NF- κ B accounts for stress-induced anxiety behaviors via enhancing nNOS-CAPON-Dexas1 coupling. *Journal of neurochemistry*, 146(5):598-612.

227. Zhu, W., Ge, X., Gao, P., Li, M., Guan, Y. and Guan, X. (2018). Adolescent cocaine exposure induces prolonged synaptic modifications in medial prefrontal cortex of adult rats. *Brain Structure and Function*, 223 (4):1829-1838.

**For more references, please visit our website at
www.fdneurotech.com.**